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hazards pictogram. Numbers below refer to performance levels

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CHLOROSULFONATED POLYETHYLENE 4/10 DRY BOX GLOVES AGAINST CHEMICAL, MICRO-ORGANISMS, MECHANICAL HAZARDS AND RADIOACTIVE CONTAMINATION.

We thank you for your confidence. Please read this brochure before using the gloves. Gloves have to be forwarded and stored in their original packing.

Gloves have to be stored in a dry place, sheltered from light, at temperatures between 5° and 25° C. Gloves should not be stored close to electric

installations. Those glove must be used in ambient temperature. Before each use , glove integrity should be checked. We remind you that the choice and

use of a Personal Protective Equipment (PPE) must correspond to the European Directive 89/656/CEE dated 30/11/89: in particular, employer must have

as a preliminary study the hazards binded to intervention and those generated by the use of an IPE (for example tightness breaking) For monting and

dismounting gloves on their port, follow the instructions written by security responsible and port manufacturer. Ensure users hands are dry, clean, with

short nails and no jewellery. There are no allergens in the glove. This glove has a good resistance to ionizing radiations but is not protecting the hands

against this hazard. If necessary, gloves can be washed with soapy water, rinsed under tap water, dried in low temperature current of air. For other

chemical products please contact us Those type of gloves can be equipped with a support ring, which can be adapted according to different diameters. In

this case, the glove's reference will be followed by the following marking: JA(ambidextrous) +BS, GJD(right) +BS, GJG(left) +BS, or AD (special bead

diameter) and A for Ambidextrous and D for right hand or G for left hand. The customer must not make installation of these support rings. Those gloves

are not concerned by external disinfection. Concerning internal disinfection, a disposable glove can be used for each handling. Life time of that glove, in

its original packing and stored in conditions recommended above, is 3 years.

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PIERCAN S.A. : 11 et 11 bis rue Charbonnel 75013 PARIS Tél : 33 (0)1 45 88 66 27 Fax : 33 (0)1 45 80 98 30